

Protocol**Immunohistochemistry (IHC)****Reagents Required**

Product	Preparation	Item No.
Phosphate Buffered Saline (PBS)	Use 10X PBS, pH 7.2 (0.2 M Potassium Phosphate, 1.5 M NaCl). Dilute appropriate volume to 1X with deionized water.	MB-008
Xylene		
95% and 100% Ethanol		
UltraPure Sterile Water		MB-009-1000
Antibody Dilution Buffer	Prepare 100 mL of PBS, supplemented with 1 mL of normal serum of same species as host for the secondary antibody.	
30% Hydrogen Peroxide Solution		KHJ001
Biotinylated Secondary Antibody, 1:500		
Streptavidin Peroxidase Conjugated, 1:500		KHD001
DAB Substrate or TMB Membrane Peroxidase Substrate for stable brown or blue staining, respectively.		DAB-10 / TMBM-100
Polymount Mounting Media		KHH001

Procedure for Frozen Sections

1. Snap-freeze fresh tissues in liquid nitrogen or isopentane pre-cooled in liquid nitrogen, embedded in OCT compound in cryomolds. Store frozen blocks at -80°C.
2. Cut cryostat sections 4–8 mm thick and mount onto Superfrost™ Plus slides or gelatin-coated slides. Store slides at -80°C until needed.
3. Before staining, warm slides at room temperature for 30 minutes and fix in ice-cold acetone for 10 minutes. Air-dry for 30 minutes.
4. Wash in PBS. Read absorbances on plate using microplate reader.

Procedure for Paraffin Sections

1. Deparaffinize sections in xylene 2 times for 5 minutes.
2. Hydrate with 100% ethanol 2 times for 3 minutes.
3. Hydrate with 95% ethanol for 1 minute.

4. Rinse in UltraPure sterile water.

Procedure for Immunoenzyme Staining

1. Follow procedure for pretreatment as required.
2. Rinse sections in PBS 2 times for 2 minutes.
3. Incubate sections in normal serum block with the same species as the secondary antibody (e.g. Normal Goat Serum (NGS) if secondary antibody is goat host).
Note: Since this protocol uses avidin-biotin detection system, avidin/biotin block may be needed based on tissue type. If you do, the avidin/biotin blocking should be done after normal serum block and before primary antibody incubation.
4. Incubate sections in primary antibody at appropriate dilution in dilution buffer for 1 hour at room temperature or overnight.
Note: Do not rinse sections between serum block and primary antibody incubation.
5. Rinse in PBS buffer 3 times for 2 minutes.
6. Incubate sections in 1% hydrogen peroxidase in PBS for 10 minutes at room temperature.
7. Rinse in PBS buffer 3 times for 2 minutes.
8. Incubate sections in biotinylated secondary antibody in PBS buffer for 30 minutes at room temperature.
9. Rinse in PBS buffer 3 times for 2 minutes.
10. Incubate sections in streptavidin peroxidase in PBS buffer for 30 minutes at room temperature.
11. Rinse in PBS buffer 3 times for 2 minutes.
12. Incubate sections in peroxidase substrate solution.
13. Rinse in PBS buffer 3 times for 2 minutes.
14. Rinse in UltraPure water 3 times for 5 minutes.
15. Dehydrate through 95% ethanol for 1 minute, 100% ethanol 2 times for 3 minutes.
16. Clear in xylene 2 times for 5 minutes.
17. Coverslip with mounting medium.

Rockland products are for research use only and are not intended for therapeutic or diagnostic applications. Please contact a technical service representative for more information. All properties listed are typical characteristics and are not specifications. All suggestions and data are offered in good faith but without guarantee as conditions and methods of use of our products are beyond our control. All claims must be made within 30 days following the date of delivery. Suggested uses of our products are not recommendations to use our products in violation of any patent or as a license under any patent of Rockland Immunochemicals, Inc. If you require a commercial license to use this material and do not have one, then return this material, unopened to: Rockland Immunochemicals, Inc., P.O. BOX 5199, Limerick, Pennsylvania 19468, USA.

© 2025 Rockland Immunochemicals, Inc. All rights reserved.